



## BIOSTAT® CultiBag RM Culturing Convenience



#2

Application  
Note

#3

Protocol for fed  
batch, serum free  
cultivation of CHO  
XM 111 suspension  
cells in the BIOSTAT  
CultiBag RM 20

#4

#5

#6

Dipl. Ing. Irina Bauer\*, Prof. Dr. Regine Eibl\*,  
Dr. Thorsten Adams\*\*

\* Zurich University of Applied Sciences  
(ZHAW), Wädenswil, Switzerland

\*\* Sartorius Stedim Biotech GmbH,  
August-Spindler-Str.11,  
37079, Goettingen, Germany

### Introduction

In this application note, we describe a protocol for the propagation of the model protein secreted alkaline phosphatase expressing *CHO XM 111* suspension cells (obtained from Prof. Dr. Martin Fussenegger, Swiss Federal Institute of Technology, Zurich) in selective, chemically defined, protein and peptide-free ChoMaster media (HP-1 and HP-5) using the disposable bioreactor BIOSTAT CultiBag RM 20 basic.

Generally, the inoculum for the bioreactor is prepared by pooling T-flasks. The pre-culture is routinely realized in 75 cm<sup>2</sup> and 175 cm<sup>2</sup> culture flasks containing CHOMaster FMX-8 growth medium, which was used for maintenance of the culture. In order to ensure optimum growth in T-flasks, the CHO suspension cells are incubated at 37°C in a humidified atmosphere of 10% CO<sub>2</sub> in air. Cells are seeded at a minimal density of 2–3 × 10<sup>5</sup> viable



Fig. 1: BIOSTAT® CultiBag RM20 basic.

cells/mL and subcultured or inoculated in the larger scale when cell densities have reached values around 1\*10<sup>6</sup> viable cells/mL.

In our experience, this method described for *CHO XM 111* suspension cells can also be successfully applied to other animal cell lines such as non-transfected CHO suspension cells, *Sf-9/Sf-21* suspension cells (DSMZ), and engineered *HEK-293 EBNA* suspension cells (Cytos Biotechnology AG, Switzerland). Modifications mainly concern the culture medium.

## 1. Equipment and Material

- BIOSTAT® CultiBag RM 20 basic (Sartorius Stedim Biotech GmbH)
- CultiBag RM 2L (with Sartofluor 300 Capsule, 0.2 µm; Sartorius Stedim Biotech GmbH)
- CHOMaster media HP-1/HP-5 (Cell Culture Technologies GmbH)
- Cedex Cell Counter (Innovatis AG) or Cedex HiRes (Innovatis AG) or NucleoCounter (ChemoMetec A/S)
- Bioprofile Analyzer 100 or BioProfile Analyzer 100+ (Nova Biomedical)
- T-flasks (T-75, T-175)
- CO<sub>2</sub> incubator
- Water bath
- Magnetic stirrer
- Pipetboy (Integra Biosciences AG)
- Peristaltic-pump (e.g. Dose-it 803, Vitaris AG)
- Sterile syringes (10mL, 50 mL)
- Serological pipettes
- Reaction tubes and sample vials (1.5 mL)
- Sterile bottles
- Sterile aluminium foil
- Safety cabinet class II
- Laminar flow module
- Roll-Boy with tripod

## 2. Methods

### a. Schedule

Day 1:	Establishment of preculture I in T-75 flask with rapidly growing, healthy <i>CHO XM 111</i> suspension cells characterized by logarithmic growth and doubling times $\leq 24$ hours
Day 3:	Feeding of preculture I with ChoMaster FMX-8 growth medium
Day 5:	Establishment of preculture II (T-175 flask) from preculture I Passage cells into T-175 (minimal seeding density of $2\text{-}3 \times 10^5$ viable cells/mL), if cell density has reached $1 \times 10^6$ viable cells/mL
Day 7:	Pooling of preculture II, inoculation and starting-up BIOSTAT CultiBag RM 20 with the disposable bioreactor bag CultiBag RM 2L operating with 100 mL cell suspension ( $1 \times 10^6$ viable cells/mL) and 100 mL ChoMaster HP-1 growth medium (see section 2d, 2e, 2f and 2g)
Day 7:	Fermentor/Bioreactor and medium preparation (see section 2b and 2c)
Day 8, 9, 10, 11:	Sampling, successive feeding of ChoMaster growth medium (up to cell densities of $1.2 \times 10^6$ viable cells/mL HP-1, subsequent feeding of HP-5 growth medium), increase of rocking rate and IPC (see section 3 and 5). The feeding procedure should be also done in such a mode that glucose levels below 1.0 g/L are avoided.
Day 12 or 13:	Partial or complete harvest of cells (see section 4). Cell densities between 2 and $4 \times 10^6$ viable cells/mL may be achievable. Aim at viabilities above 95%.

### b. Fermentor|Bioreactor preparation

100 mL of ChoMaster HP-1 growth medium containing 0.2 % Pluronic are filled in the CultiBag RM 2L in the safety cabinet (clamped air filters)

*Keep in mind that there is no need to use Pluronic in media containing serum.*

### c. Media

Selective medium for T-flasks: filter-sterilized, conditioned (37 °C, pH 7.3) ChoMaster FMX-8 growth medium (Cell Culture Technologies).

Additional supplements for FMX-8 medium:

Used antibiotics to keep cells under selection pressure, support cell growth and prevent SEAP expression: 100 mg mL<sup>-1</sup> G418 sulphate, 5 mg mL<sup>-1</sup> puromycin dihydrochloride, 2.5 mg mL<sup>-1</sup> tetracycline hydrochloride.

Medium for BIOSTAT CultiBag RM 20: filter-sterilized, conditioned (37 °C, pH 7.3) ChoMaster HP-1- and HP-5 growth medium (Cell Culture Technologies)

Additional supplements for HP-1 and HP-5 medium:

2.5 mg mL<sup>-1</sup> tetracycline hydrochloride, supports cell growth and prevents SEAP expression and 0.2 % Pluronic F68 solution (Sigma) protects cells against shear (*only necessary in serum-free media!*)

### d) Preculture and Inoculum for CultiBag RM 2L

For establishing the preculture II (T-175) representing the subsequent inoculum after pooling procedure approximately 4 days are required. In case of use of cryopreserved vials we recommend a previous T-flask cultivation of 14 days. In other words, the use of cryopreserved vials instead of T-75 will prolong the precultivation time.

In order to obtain the desired seeding cell density of about 5\*10<sup>5</sup> viable cells/mL for the CultiBag RM 2L, harvest of 5\*10<sup>7</sup> viable cells from T-flasks, pooling of the cell pellets and resuspension in 100 mL fresh ChoMaster HP-1 growth medium have to be carried out.

Consequently, the cells were transferred from T-175 into a sterile beaker (pipetting) covered with a sterile aluminum foil and

incubated (CO<sub>2</sub> incubator) for 3 hours in order to allow the cells to settle. Alternatively for other cell lines, the cells can be centrifuged at maximum 200 g.

The consumed growth medium (FMX-8) was then aspirated and replaced with fresh HP-1 growth medium (pH 7.3, 37°C) in the safety cabinet. After cell density check the cell suspension in the sterile beaker was ready for its use in CultiBag RM 2L.

### e. Corrective agent

Acid:	0-10% CO <sub>2</sub>
-------	-----------------------

### f. Culture conditions

Starting culture volume:	200 mL
Final culture volume:	1000 mL
Rocking rate:	14-30 rpm
Rocking angle:	6°
pH:	7.3-6.9
Temperature:	37 °C
Aeration rate:	0.2 vvm
Start cell density:	5 x 10 <sup>5</sup> viable cells/mL
Final cell density:	3-4 x 10 <sup>6</sup> viable cells/mL
Cultivation time:	6-7 days

### g. Inoculation

- By inserting a syringe into the CultiBag's luer lock inoculation port, 100 mL of the prepared cell suspension (see section 2d) were added in the safety cabinet (exhaust air filter was clamped off).
- The filled CultiBag (100 mL HP-1 growth medium, see section 2b, and 100 mL cell suspension, see section 2d) was put back on the tray (clamped air filters), fixed and installed.
- The filter heater was installed and switched on.
- Air filter lines were opened and aeration (0.2 vvm), rocking (14 rpm, 6°) and heating (37 °C) were switched on.

### 4. Complete cell harvest or Scale-up

For harvesting the cells/product, one of the attached ports can be used. For the scale-up into a larger volume the following procedure can be used:

- The BIOSTAT CultiBag RM basic station was switched off.
- The air filters were closed.
- The CultiBag RM 2L was removed from the tray and transferred to a laminar flow module.
- The CultiBag was hung on a tripod standing on a Roll-Boy in the laminar flow module.
- The exhaust filter of the CultiBag was opened whereas inlet filter was closed.

### 3. Start-up and operation of BIOSTAT CultiBag RM 20

½ hours:

Sample 0: Analytics (section 5)

---

24 hours:

Sample 1: Analytics (section 5)

---

48 hours:

Sample 2: Analytics (section 5), feeding with 200 mL HP-1 growth medium and rocking rate increase (16 rpm)

---

72 hours:

Sample 3: Analytics (section 5), feeding with 200 mL HP-5 and rocking rate increase (18 rpm)

---

96 hours:

Sample 4: Analytics (section 5), feeding with 200 mL HP-5 and rocking rate increase 25 rpm

---

120 hours:

Sample 5: Analytics (section 5), feeding with 200 mL HP-5 and rocking rate increase 30 rpm

---

6 or 7 days:

Sample 6 and 7: Analytics (section 5), partial or complete cell harvest (section 4). In case of partial cell suspension harvest, the adequate amount of fresh HP-5 growth medium is fed.

- 200 mL of HP-5 growth medium were fed.

- For about 3 hours the cells were allowed to settle on the bottom of the CultiBag RM 2L.

- The medium was removed from the CultiBag using the tube of the fill/harvest port.

### 5. Analytics

Daily one 2 mL sample is taken in order to determine:  
Cell growth and viability (1 mL sample) by use of Cedex or NucleoCounter instead of traditional, time-consuming, manual cell counting (hemocytometer, Trypan Blue)  
Glucose, lactate, glutamine, glutamate, pH (1 mL sample) by use of Nova BioProfile Analyzer 100 or its successor. Alternatively, other automatized analyzers (e.g. YSI 2700 Bio-chemistry Analyzer, YSI Incorporated, and Eppendorf Ebio plus) or also test kits (for example, from Roche Diagnostics) are available.

# Sales and Service Contacts

For further contacts, visit [www.sartorius-stedim.com](http://www.sartorius-stedim.com)

## Europe

### Germany

Sartorius Stedim Biotech GmbH  
August-Spindler-Strasse 11  
37079 Goettingen

Phone +49.551.308.0  
Fax +49.551.308.3289

[www.sartorius-stedim.com](http://www.sartorius-stedim.com)

Sartorius Stedim Systems GmbH  
Schwarzenberger Weg 73-79  
34212 Melsungen

Phone +49.5661.71.3400  
Fax +49.5661.71.3702

[www.sartorius-stedim.com](http://www.sartorius-stedim.com)

### France

Sartorius Stedim Biotech S.A.  
Z.I. des Paluds  
Avenue de Jouques – BP 1051  
13781 Aubagne Cedex

Phone +33.442.845600  
Fax +33.442.845619

Sartorius Stedim France SAS  
ZI des Paluds  
Avenue de Jouques – CS 71058  
13781 Aubagne Cedex

Phone +33.442.845600  
Fax +33.442.846545

### Austria

Sartorius Stedim Austria GmbH  
Franzosengraben 12  
A-1030 Vienna

Phone +43.1.7965763.18  
Fax +43.1.796576344

### Belgium

Sartorius Stedim Belgium N.V.  
Leuvensesteenweg, 248/B  
1800 Vilvoorde

Phone +32.2.756.06.80  
Fax +32.2.756.06.81

### Denmark

Sartorius Stedim Nordic A/S  
Hoerskaetten 6D, 1.  
DK-2630 Taastrup

Phone +45.7023.4400  
Fax +45.4630.4030

### Italy

Sartorius Stedim Italy S.p.A.  
Via dell'Antella, 76/A  
50012 Antella-Bagno a Ripoli (FI)

Phone +39.055.63.40.41  
Fax +39.055.63.40.526

### Netherlands

Sartorius Stedim Netherlands B.V.  
Edisonbaan 24  
3439 MN Nieuwegein

Phone +31.30.6025080  
Fax +31.30.6025099

### Spain

Sartorius Stedim Spain SA  
C/Isabel Colbrand 10-12,  
Planta 4, Oficina 121  
Polígono Industrial de Fuencarral  
28050 Madrid

Phone +34.91.3586102  
Fax +34.91.3588804

### Switzerland

Sartorius Stedim Switzerland GmbH  
Lerzenstrasse 21  
8953 Dietikon

Phone +41.44.741.05.00  
Fax +41.44.741.05.09

### U.K.

Sartorius Stedim UK Limited  
Longmead Business Park  
Blenheim Road, Epsom  
Surrey KT19 9 QQ

Phone +44.1372.737159  
Fax +44.1372.726171

## America

### USA

Sartorius Stedim North America Inc.  
5 Orville Drive  
Bohemia, NY 11716

Toll-Free +1.800.368.7178  
Fax +1.631.254.4253

Sartorius Stedim SUS Inc.  
1910 Mark Court  
Concord, CA 94520

Phone +1.925.689.6650  
Toll Free +1.800.914.6644  
Fax +1.925.689.6988

Sartorius Stedim Systems Inc.  
201 South Ingram Mill Road  
Springfield, MO 65802

Phone +1.417.873.9636  
Fax +1.417.873.9275

### Argentina

Sartorius Argentina S.A.  
Int. A. Avalos 4251  
B1605ECS Munro  
Buenos Aires

Phone +54.11.4721.0505  
Fax +54.11.4762.2333

### Brazil

Sartorius do Brasil Ltda  
Av. Dom Pedro I, 241  
Bairro Vila Pires  
Santo André  
São Paulo  
Cep 09110-001

Phone +55.11.4451.6226  
Fax +55.11.4451.4369

### Mexico

Sartorius de México S.A. de C.V.  
Circuito Circunvalación Poniente No. 149  
Ciudad Satélite  
53100 Naucalpan, Estado de México

Phone +52.5555.62.1102  
Fax +52.5555.62.2942

## Asia | Pacific

### China

Sartorius Stedim Beijing  
Representative Office  
No. 33, Yu'an Road,  
Airport Industrial Zone B, Shunyi District  
Beijing 101300

Phone +86.10.80426516  
Fax +86.10.80426580

Sartorius Stedim Shanghai  
Representative Office  
Room 618, Tower 1, German Centre,  
Shanghai, PRC., 201203

Phone +86.21.28986393  
Fax +86.21.28986392.11

Sartorius Stedim Guangzhou Office  
Room 704, Broadway Plaza,  
No. 233-234 Dong Feng West Road  
Guangzhou 510180

Phone +86.20.8351.7921  
Fax +86.20.8351.7931

### India

Sartorius Stedim India Pvt. Ltd.  
10, 6th Main, 3rd Phase Peenya  
KIADB Industrial Area  
Bangalore – 560 058

Phone +91.80.2839.1963|0461  
Fax +91.80.2839.8262

### Japan

Sartorius Stedim Japan K.K.  
KY Building, 8-11  
Kita Shinagawa 1-chome  
Shinagawa-ku  
Tokyo 140-0001

Phone +81.3.3740.5407  
Fax +81.3.3740.5406

### Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.  
Lot L3-E-3B, Enterprise 4  
Technology Park Malaysia  
Bukit Jalil  
57000 Kuala Lumpur

Phone +60.3.8996.0622  
Fax +60.3.8996.0755

### Singapore

Sartorius Stedim Singapore Pte. Ltd.  
10, Science Park Road, The Alpha  
#02-25, Singapore Science Park 2  
Singapore 117684

Phone +65.6872.3966  
Fax +65.6778.2494

### Australia

Sartorius Stedim Australia Pty. Ltd.  
Unit 5, 7-11 Rodeo Drive  
Dandenong South Vic 3175

Phone +61.3.8762.1800  
Fax +61.3.8762.1828